

KERN & Sohn GmbH

Ziegelei 1 D-72336 Balingen E-mail:info@kern-sohn.com Tel: +49-[0]7433- 9933-0 Fax: +49-[0]7433-9933-149 Internet: www.kern-sohn.com

Operating Manual Precision balance



Version 1.5 02/2013 GB



PCB-BA-e-1315



KERN PCB

Version 1.5 02/2013 Operating Manual Compact balance

Table of Contents

1	Technical Data	4
2	Declaration of conformity	7
3 3.1 3.2 3.3 3.4	Basic Information (General) Proper use Improper Use Warranty Monitoring of Test Resources	8 8 8
4 4.1 4.2	Basic Safety Precautions Pay attention to the instructions in the Operation Manual Personnel training	9
5 5.1 5.2	Transportation & Storage Testing upon acceptance Packaging	9
6 6.1 6.2 6.3 6.4 6.5 6.6 6.7 6.8	Unpacking, Setup and Commissioning Installation Site, Location of Use Unpacking/erection Mains connection Operation using a (rechargeable) battery (optional) Connection of peripheral devices Initial Commissioning Adjustment	9 10 11 11 12
6.9	Underfloor weighing	13
7 7.1 7.2	Operation elements	14
	Basic Operation	15
Taring. PRE-T	ng ARE-Function	5 6
Parts c	inus weighings ounting al weighings	17
Percen	t determination	19

9	Menu	
9.1	Navigation in the menu	
9.2	Menu overview	
9.3	Description of individual menu items	
Weigh	ing Units	
Dosing	g and Zero-tracking	
Select	ion adjustment weight	
	y background illumination	
Anima	I weighing function	
Reset	to factory setting	
9.4	Interface parameters	
Data tr	ransfer mode	
printo	ut	
Baud r	rate	
10	Data output RS 232 C	
10.1	Technical Data	
10.2	Pin allocation of balance output bushing:	
10.3	Explanation of the data transfer	
10.4	Output on bar code printer	
11	Service, maintenance, disposal	
11.1	CLEANING	
11.2	Service, maintenance	
11.3	Disposal	
12	Instant help	

1 Technical Data

KERN	PCB 100-3	PCB 200-2	PCB 250-3	PCB 350-3
Readability (d)	0,001 g	0,01 g	0,001 g	0,001 g
Weighing range (max) (d)	100 g	200 g	250 g	350 g
Taring range (subtractive)	100 g	200 g	250 g	350 g
Reproducibility	0,001 g	0,01 g	0,001 g	0,002 g
Linearity	±0,003 g	± 0,02 g	±0,003 g	±0,004 g
Minimum unit weight at piece counting	0,002 g	0,02 g	0,002 g	0,002 g
Warm-up time		2 ho	ours	
Reference quantities at piece counting		5, 10, 20	, 25, 50	
Weighing Units	Detail	s " Weighing ι	inits " see ch	ot. 9.3
Recommended adjustment weight, not added (class) Details for "Selection of the Adjustment weight" in ch. 9.3	100 g (F1)	200 g (M1)	200 g (F1)	300 g (F1)
Stabilization time (typical)	3 sec.			
Operating temperature	+ 5° C + 35° C			
Humidity of air		max. 80 % (no	t condensing)	
Housing (B x D x H) mm		163 x 24	45 x 79	
Windscreen mm	Ø 106, high 42 - Ø 106, high 42		high 42	
Weighing plate mm	Ø 81	Ø 105	Ø 105 Ø 81	
Weight kg (net)	1,1			
Electric Supply	220V-240V AC, 50 Hz / 9 V, 300 mA			
Battery operation	9 V- compound battery (optional)			
Rechargeable battery (optional)	Operating time 24 h with display background illumination Operating time 48 h without display background illumination/loading time 8 h			
Interface	RS 232			
Underfloor weighing	standard			

KERN	PCB 1000-1	PCB 1000-2	PCB 2000-1	PCB 2500-2	
Readability (d)	0,1 g	0,01 g	0,1 g	0,01 g	
Weighing range (max) (d)	1000 g	1000 g	2000 g	2500 g	
Taring range (subtractive)	1000 g	1000 g	2000 g	2500 g	
Reproducibility	0,1 g	0,01 g	0,1 g	0,01 g	
Linearity	±0,2 g	±0,03 g	±0,2 g	±0,03 g	
Minimum unit weight at piece counting	0,2 g	0,02 g	0,2 g	0,02 g	
Warm-up time	30 minutes	2 hours	30 minutes	2 hours	
Reference quantities at piece counting		5, 10, 20), 25, 50		
Weighing Units	Detai	ls " Weighing	units" see chp	ot. 9.3	
Recommended adjustment weight, not added (class) Details for "Selection of the Adjustment weight" in ch. 9.3	1000 g (M1)	1000 g (F1)	2000 g (M1)	2000 g (F1)	
Stabilization time (typical)	3 sec.				
Operating temperature	+ 5° C + 35° C				
Humidity of air		max. 80 % (no	ot condensing)		
Housing (B x D x H) mm		163 x 2	45 x 79		
Weighing plate mm	130 x 130				
Weight kg (net)	1,4				
Electric Supply	220	V-240V AC, 50) Hz / 9 V, 300	mA	
Battery operation	9	V- compound I	pattery (optiona	al)	
Rechargeable battery (optional)	Operating time 24 h with display background illumination Operating time 48 h without display background illumination/loading time 8 h				
Interface	RS 232				
Underfloor weighing	standard				

KERN	PCB 3500-2	PCB 6000-1	PCB 6000-0	PCB 10000-1
Readability (d)	0,01 g	0,1 g	1 g	0,1 g
Weighing range (max) (d)	3500 g	6 kg	6 kg	10 kg
Taring range (subtractive)	3500 g	6 kg	6 kg	10 kg
Reproducibility	0,02 g	0,1 g	1 g	0,1 g
Linearity	0,04 g	± 0,3 g	± 2 g	± 0,3 g
Minimum unit weight at piece counting	0,02 g	0,2 g	2 g	0,2 g
Warm-up time	2 hours	2 hours	30 minutes	2 hours
Reference quantities at piece counting		5, 10, 20), 25, 50	
Weighing Units	Details "Weighing units" see chpt. 9.3			t. 9.3
Recommended adjustment weight, not added (class) Details for "Selection of the Adjustment weight" in ch. 9.3	3000 g (F1)	5000 g (F2) 5000 g (M		10 kg (F1)
Stabilization time (typical)		3 s	ec.	
Operating temperature		+ 5° C	. + 35° C	
Humidity of air		max. 80 % (no	ot condensing)	
Housing (B x D x H) mm		163 x 2	45 x 79	
Weighing plate mm	130 x 130 150 x 170 150 x		150 x 170	150 x 170
Weight kg (net)	1.4	2.0	2.0	2.0
Electric Supply	220	V-240V AC, 50) Hz / 9 V, 300	mA
Battery operation	9 V- compound battery (optional)			al)
Rechargeable battery (optional)	Operating time 24 h with display background illumination Operating time 48 h without display background illumination/loading time 8 h			
Interface	RS 232			
Underfloor weighing	standard			

2 Declaration of conformity



KERN & Sohn GmbH D-72322 Balingen-Frommern Postbox 4052 email: info@kern-sohn.de

Phone: 0049-[0]7433- 9933-0 Fax: 0049-[0]7433-9933-149 Internet: www.kern-sohn.de

Declaration of conformity

EC- D EC-Di EC- D	onformitätserklä Déclaration de co ichiarazione di co Declaração de co eklaracja zgodn	onformité EC-Declaración de Conformidad conformità EC-Conformiteitverklaring onformidade EC- Prohlášení o shode		
D	Konformitäts- erklärung	Wir erklären hiermit, dass das Produkt, auf das sich diese Erklärung bezieht, mit den nachstehenden Normen übereinstimmt.		
GB	Declaration of conformity	We hereby declare that the product to which this declaration refers conforms to the following standards.		
CZ	Prohlášení o shode	Tímto prohlašujeme, že výrobek, kterého se toto prohlášení týká, je v souladu s níže uvedenými normami.		
Е	Declaración de conformidad	Manifestamos en la presente que el producto al que se refiere esta declaración está de acuerdo con las normas siguientes		
F	Déclaration de conformité	Nous déclarons avec cela responsabilité que le produit, auquel se rapporte la présente déclaration, est conforme aux normes citées ci-après.		
I	Dichiarazione di conformitá	Dichiariamo con ciò che il prodotto al quale la presente dichiarazione si riferisce è conforme alle norme di seguito citate.		
NL	Conformiteit- verklaring	Wij verklaren hiermede dat het product, waarop deze verklaring betrekking heeft, met de hierna vermelde normen overeenstemt.		
Ρ	Declaração de conformidade	Declaramos por meio da presente que o produto no qual se refere esta declaração, corresponde às normas seguintes.		
PL	Deklaracja zgodności	Niniejszym oświadczamy, że produkt, którego niniejsze oświadczenie dotyczy, jest zgodny z poniższymi normami.		
RUS	Заявление о соответствии	Мы заявляем, что продукт, к которому относится данная декларация, соответствует перечисленным ниже нормам.		

Electronic Balance: KERN FCB, FCE, FKB...A, PCB

EU Directive	Standards
2004/108/EC	EN 61326-1: 2006
	EN 61326-2-2: 2006
	EN 61000-3-2: 2006
	EN 61000-3-3: 2008
2006/95/EC	EN 60950-1:2006+A11:2009

Suntin

Signature: ______KER

KERN & Sohn GmbH Management

KERN & Sohn GmbH, Ziegelei 1, D-72336 Balingen, Tel. +49-[0]7433/9933-0, Fax +49-[0]7433/9933-149

Date: 06.03.2012

3 Basic Information (General)

3.1 Proper use

The balance you purchased is intended to determine the weighing value of material to be weighed. It is intended to be used as a "non-automatic" balance, i.e. the material to be weighed is manually and carefully placed in the centre of the weighing plate. As soon as a stable weighing value is reached the weighing value can be read.

3.2 Improper Use

Do not use balance for dynamic weighing. In the event that small quantities are removed or added to the material to be weighed, incorrect weighing results can be displayed due to the "stability compensation" in the balance. (Example: Slowly draining fluids from a container on the balance). Do not leave permanent load on the weighing plate. This may damage the measuring system. Impacts and overloading exceeding the stated maximum load (max) of the balance, minus a possibly existing tare load, must be strictly avoided. Balance may be damage by this.

Never operate balance in explosive environment. The serial version is not explosion protected.

The structure of the balance may not be modified. This may lead to incorrect weighing results, safety-related faults and destruction of the balance.

The balance may only be used according to the described conditions. Other areas of use must be released by KERN in writing.

3.3 Warranty

Warranty claims shall be voided in case

- Our conditions in the operation manual are ignored
- The appliance is used outside the described uses
- The appliance is modified or opened
- Mechanical damage and damage caused by media, liquids
- Natural wear and tear
- The appliance is improperly set up or incorrectly electrically connected
- The measuring system is overloaded

3.4 Monitoring of Test Resources

In the framework of quality assurance the measuring-related properties of the balance and, if applicable, the testing weight, must be checked regularly. The responsible user must define a suitable interval as well as type and scope of this test. Information is available on KERN's home page (www.kern-sohn.com) with regard to the monitoring of balance test substances and the test weights required for this. In KERN's accredited DKD calibration laboratory test weights and balances may be calibrated (return to the national standard) fast and at moderate cost.

4 Basic Safety Precautions

4.1 Pay attention to the instructions in the Operation Manual



Carefully read this operation manual before setup and commissioning, even if you are already familiar with KERN balances.

Versions in other languages are non-binding translations. The only binding version is the original document in German.

4.2 Personnel training

The appliance may only be operated and maintained by trained personnel.

5 Transportation & Storage

5.1 Testing upon acceptance

When receiving the appliance, please check packaging immediately, and the appliance itself when unpacking for possible visible damage.

5.2 Packaging

Keep all parts of the original packaging in case you need to return the appliance. Only use original packaging for returning.

Before sending, disconnect all connected cables and loose/movable parts. Attach possibly existing transport safeguards. Secure all parts, e.g. weighing plate, mains adapter etc., to prevent slipping and damage.

6 Unpacking, Setup and Commissioning

6.1 Installation Site, Location of Use

The balances are designed in a way that reliable weighing results are achieved in common conditions of use.

You will work accurately and fast, if you select the right location for your balance. **Therefore, observe the following for the installation site:**


- Place the balance on a firm, level surface;
- Avoid extreme heat as well as temperature fluctuation caused by installing next to a radiator or in the direct sunlight;
- Protect the balance against direct draughts due to open windows and doors;
- Avoid jarring during weighing;
- Protect the balance against high humidity, vapors and dust;
- Do not expose the device to extreme dampness for longer periods of time. Nonpermitted condensation (condensation of air humidity on the appliance) may occur if a cold appliance is taken to a considerably warmer environment. In this case, acclimatize the disconnected appliance for ca. 2 hours at room temperature.
- Avoid static charge of goods to be weighed and weighing container.

Major display deviations (incorrect weighing results) may be experienced should electromagnetic fields (e.g. due to mobile phones or radio equipment), static electricity accumulations or instable power supply occur. Change location or remove source of interference.

6.2 Unpacking/erection

Carefully remove the balance from the packaging, remove plastic cover and setup balance at the intended workstation.

The balance must be installed in a way that the weighing plate is exactly in horizontal position.

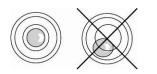

Level balance with foot screws until the air bubble of the water balance is in the prescribed circle.

Scope of delivery / serial accessories

- Balance
- Weighing plate
- Mains power supply
- Protective cover
- Operating Manual
- Windscreen (PCB 100-3, PCB 250-3, PCB 350-3)

6.3 Mains connection

Power is supplied via the external mains adapter. The stated voltage value must be the same as the local voltage.

Only use original KERN mains adapters. Using other makes requires consent by KERN.

6.4 Operation using a (rechargeable) battery (optional)

Lift-off the battery cover on the lower side of the balance. Connect 9 V compound battery.

Replace the battery compartment cover.

For battery operation the balance has an automatic switch-off function which can be activated or deactivated in the menu (chapter 9).

- ⇒ In weighing mode keep the **PRINT**-key pressed until [Unit] appears.
- \Rightarrow Press **MODE** key repeatedly until "AF" appears.
- \Rightarrow Use the **SET** key, to confirm.
- \Rightarrow Use the **MODE** key to choose between the two following settings:
 - **"AF on":** In order to save the battery, the balance switches automatically off after 3 minutes without weighing.
 - "AF off": Switch-off function deactivated.
- \Rightarrow Use the **SET** key to confirm. The balance returns to weighing mode.

If the batteries are run down, "LO" appears in the display. Press **ON/OFF**-key and replace the batteries immediately.

If the balance is not used for a longer time, take out the batteries and store them separately. Leaking battery liquid could damage the balance.

If there exists an optional rechargeable battery, it has to be connected in the battery compartment via a separate plug-in socket. Now the mains adapter delivered with the rechargeable battery must be applied.

6.5 Connection of peripheral devices

Before connecting or disconnecting of additional devices (printer, PC) to the data interface, always disconnect the balance from the power supply. With your balance, only use accessories and peripheral devices by KERN, as they are ideally tuned to your balance.

6.6 Initial Commissioning

In order to obtain exact results with the electronic balances, your balance must have reached the operating temperature (see warming up time chap. 1). During this warming up time the balance must be connected to the power supply (mains, accumulator or battery).

The accuracy of the balance depends on the local acceleration of gravity. Strictly observe hints in chapter Adjustment.

6.7 Adjustment

As the acceleration value due to gravity is not the same at every location on earth, each balance must be coordinated - in compliance with the underlying physical weighing principle - to the existing acceleration due to gravity at its place of location (only if the balance has not already been adjusted to the location in the factory). This adjustment process must be carried out for the first commissioning, after each change of location as well as in case of fluctuating environment temperature. To receive accurate measuring values it is also recommended to adjust the balance periodically in weighing operation.

6.8 Adjustment

The adjustment should be made with the recommended adjustment weight (see chap. 1 "Technical data"). Adjustment is also possible with the weights of other nominal values (see table 1), but not the optimum for measuring technique.

Procedure when adjusting:

Observe stable environmental conditions. A warming up time (see chapter 1) is required for stabilization.

- \Rightarrow Turn on balance by pressing the **ON/OFF** key.
- Press the MODE key and keep it pressed, in the display appears shortly "CAL". After that the exact size appears flashing in the display (see chapter.9.3) of the adjustment weight.
- \Rightarrow Now set the adjusting weight in the centre of the weighing plate.
- ⇒ Press the SET key. Short time later there appears "CAL F", then the automatic return to the weighing mode. In the display there appears the value of the adjustment weight.

An error during adjustment or the use of an incorrect adjusting weight will result in an error message **"CAL E**". Repeat adjustment.

Keep the adjustment close to the balance. Daily control of the weighing exactness is recommended for quality-relevant applications.

6.9 Underfloor weighing

Objects unsuitable for placing on the weighing scale due to size or shape may be weighed with the help of the flush-mounted platform.

Proceed as follows:

- Switch off the balance
- Open closing cover at the balance bottom.
- Suspend hook for underfloor weighing carefully and completely.
- Place weighing balance over an opening.
- Attach load to hook and carry out weighing procedure.



Fig. 1: Setup of balance for underfloor weighing




- Always ensure that all suspended objects are stable enough to hold the desired goods to be weighed safely (danger of breaking).
- Never suspend loads that exceed the stated maximum load (max) (danger of breaking)
- Always ensure that there are no persons, animals or objects that might be damaged underneath the load.



After completing the underfloor weighing the opening on the bottom of the balance must always be closed (dust protection).

7 Operation elements

7.1 Overview of display

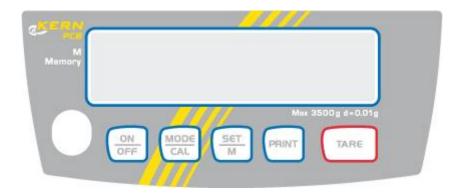

7.2 Keyboard overview

Кеу	Designation	Function		
		Calculate weighing data via interface		
PRINT	PRINT-key	Call up menu (keep key pressed until UNIT appears)		
SET	CET kov	Confirm settings in the menu		
	SET key	Save and exit menu		
		How to select menu items		
	MODE key	Change settings in the menu		
_		Adjustment		
TARE	TARE key	• Taring		
	ON/OFF switch	Turn on/off		


8 Basic Operation

Start-up




⇒ Press **ON** button.

The balance will carry out a self-test The balance is ready for weighing when the weight display appears.

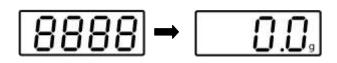

Switching Off



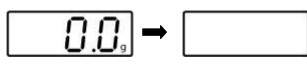

⇒ Press **OFF** button, the display disappears

Weighing

- \Rightarrow Place goods to be weighed on balance
- Wait for standstill control, after the standstill control, the weighing unit appears right hand in the display (e.g. g or kg)
- \Rightarrow Read weighing result.

If the goods are heavier than the weighing range, the display will show "**Error**" (=Overload), and a whistle is sounded.

Taring

⇒ Place an empty weighing container, the weight of the weighing container will be displayed.






⇒ Press the TARE button, the zero display disappears. The tare weight is saved until it is deleted.

00
U.U,

 \Rightarrow Weigh the material, the net weight will be indicated.

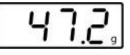

The taring process can be repeated any number of times, e.g. when adding several components for a mixture (adding). The limit is reached when the whole weighing range is exhausted.

The weight of the weighing container will be displayed as a minus number after removing the weighing container.

The tare weight is saved until it is deleted.

- C C
- ⇒ Unload the balance and press the TARE button, the zero display appears.

00
 9

PRE-TARE-Function

Delete tare

TARE



Delete PRE-TARE value



Using this function the weight of a tare vessel is stored. Even after turning off/on the weighing balance will continue working with the saved tare value.

- ⇒ In weighing mode put tare vessel on the weighing plate
- ⇒ Press repeatedly the MODE key until "PtArE" flashing appears.
- ⇒ Use SET key to store the current weight on the weighing plate as a PRE-TARE value.
- ⇒ Remove all loads from the balance and press repeatedly the MODE key until "PtArE" flashing appears.
- ⇒ Use the SET key to confirm. The PRE-TARE value is deleted, the zero display appears.

Plus/minus weighings



For example unit weight control, fabrication control etc.

- ⇒ Put the nominal weight on the weighing plate and tare using the TARE button.
- ⇒ Remove the nominal weight
- ⇒ Put the test objects subsequently on the weighing plate, the respective deviation from the nominal weight is displayed with the respective sign to "+" and "-".

According to the same procedure also packages with the same weight can be produced, referring to a nominal weight.

⇒ Back to weighing mode by pressing the **TARE** button.

Parts counting During piece counting parts can either be counted into a container or out of a container. To count a greater number of parts the average weight per part has to be determined with a small quantity (reference quantity).

The larger the reference quantity, the higher the counting exactness. High reference must be selected for small parts or parts with considerably different sizes.

The larger the reference quantity, the more accurate the parts counting.

The process has four steps:

Tare the weighing container Determine the reference unit Original weighing of reference weight Count the items



- ⇒ In weighing mode press MODE key shortly. Reference piece number "5^{PCS}" appears flashing.
- ⇒ By pressing the MODE button several times other reference quantities 5, 10, 20, 25 and 50 can be called up. Place as many pieces to count on the weighing plate as the set reference quantity requires.
- ⇒ Use the **SET** key to confirm. The balance is now in parts counting mode counting all units on the weighing plate.

Back to weighing mode by pressing the MODE button.
Error message "Er 1" Piece below minimum weight of piece (See chpt. 1 "Technical specifications"): Press MODE key and restart reference determination.
Taring The tare vessels can also be used for piece counting. Before starting piece counting use the TARE button to tare out the container.

It is useful if a mixture of several components is weighed into a tare vessel and finally the sum weight of all weighed components is necessary for control purposes (net-total, i.e. the weight of the tare vessel).

Example:

- 1. Place tare container on the weighing plate. Press the **TARE** button, the zero display disappears.
- Weigh-in component ●. Press the SET button, the zero display disappears. [▲] is displayed on the right border of the display.
- 3. Weigh-in component **②** and press **SET** key. Net-total (sum weight of the components **①** and **②**) is displayed.
- 4. Press the **SET** button, the zero display disappears.
- 5. Weigh-in component € and press SET key. Net-total (sum weight of the components € and € and €.) is displayed.
- If necessary, also fill the formula up to the desired final value. For every component more repeat the steps 4-5.
- \Rightarrow Back to weighing mode by pressing the **TARE** button.

Net-total

weighings

SET

Μ

Percent determination






Percentage calculation facilitates weight display in percent related to a reference weight equivalent to 100 %.


- ⇒ In weighing mode press MODE key repeatedly, until [100 %] is displayed.
- \Rightarrow Put a reference weight which corresponds to 100 %.
- ⇒ Store by pressing the **SET** key. Remove reference weight.
- Place goods to be weighed on balance. The weight of the sample is displayed in percentage in terms of the reference weight.


Back to weighing mode by pressing the **MODE** button.

9 Menu

9.1 Navigation in the menu

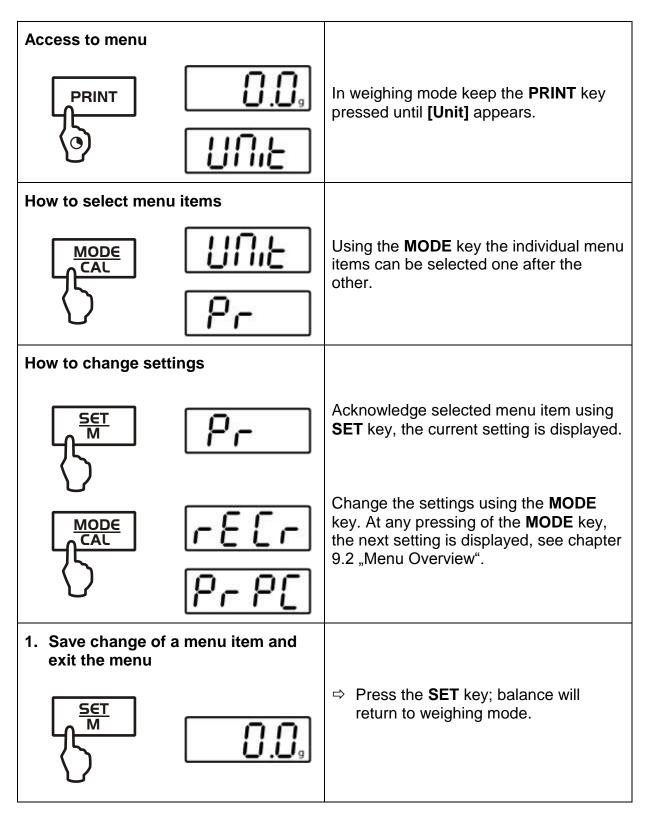

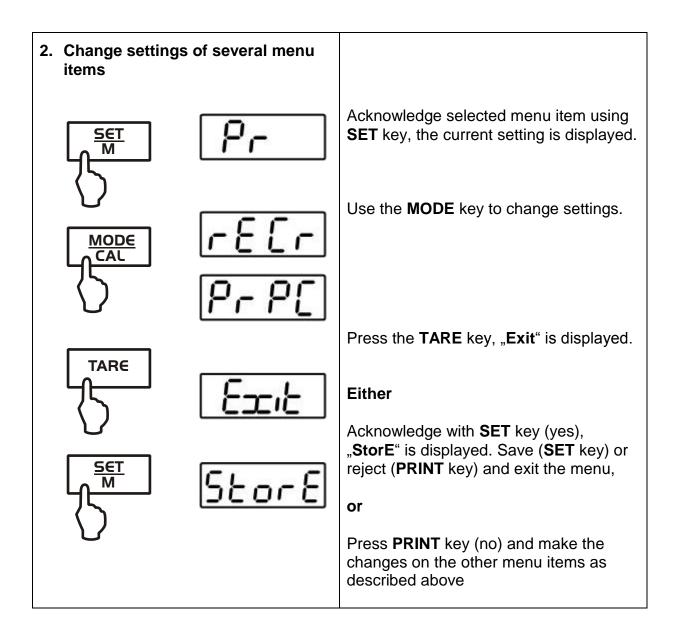

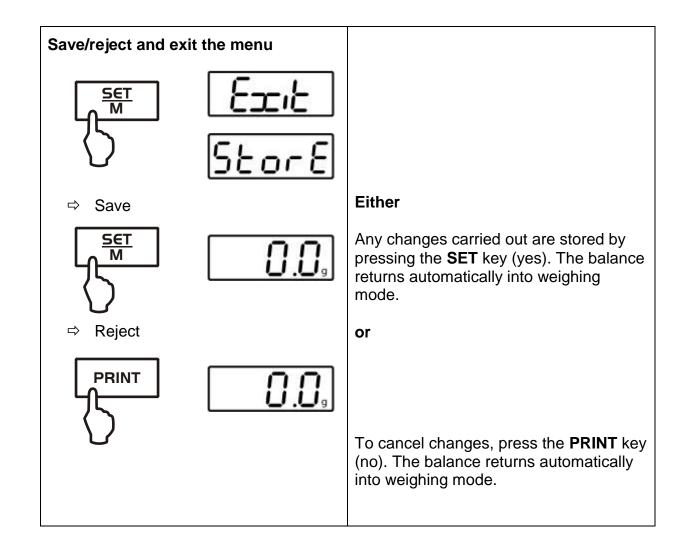

9.2 Menu overview

Description of function	Function	Parameter	Description of options
Weighing units switching	UNIT	g*	Gram
over (see chapter 9.3).		kg	Kilogram (dependent on model)
		oz	Pound
		ozt	Ounce
		lb	Troy ounce
		tlh	Tael Hongkong
		tlt	Tael Taiwan
		gn	Grain (dependent on model)
		dwt	Pennyweight (dependent on model)
		mo	Momme
		Tol	Tola
		ct	Carat (dependent on model)
		FFA	Freely selectable factor
Data transfer mode (see	Pr	rE CR*	Data output via remote control
chapter 9.4)			commands (see chapter 10.3) Data output by pressing the
		Pr PC	PRINT key (see chapter 10.3)
		AU PC	Continuous data output (see chapter 10.3)
		bA Pr	Printout on barcode printer (see chapter 10.4)
		AU Pr	Autom. Data output of stable weighing values (see chapter 10.3)
Selection printer output	LAPr	Hdr*	Edition of the headlines
(see chapter 9.4)		GrS Net	Edition of the total weight Edition of the net weight
		tAr	Edition of the tare weight
		N7E	Edition of the stored weight
		PCS	Edition of quantity
		AUJ	Edition of the unit weight
		Rqt	Edition of the reference quantity Edition of a page feeding
		FFd	at start printer output
		FFE	Edition of a page feeding
		FFE	at end printer output


Baud rate	bAUd	19200	
(see chapter 9.4).		9600*	
		4800	
		2400	
		1200	
Auto off (battery operation), see chap. 6.4	AF	on*	Automatic switch-off function after 3 min without changing load ON
		off	Automatic switch-off function after 3 min without changing load OFF
Auto Zero	tr	on*	On
(see chapter 9.3)		off	Off
Selection adjustment	CAL	100	
weight (see chapter 9.3)		200	*dependent on model
		300	
Background illumination	bL	on*	Background illumination on
of the display, (see chap. 9.3)		off	Background illumination off
0.0)		СН	The background illumination will be switched off automatically 10 sec after having reached a stable weighing value.
Animal weighing function	ANL	off*	Off
(see chapter 9.3)		3	Period 3 seconds
		5	Period 5 seconds
		10	Period 10 seconds
		15	Period 15 seconds
Reset to factory setting	rSt	no*	no
(see chap. 9.3)		yes	yes

* = default setting

9.3 Description of individual menu items

Weighing Units

⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.

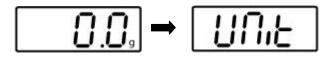

- ⇒ Press **SET** key, the selected unit is displayed.
- ⇒ Use the MODE key to select between the different units (see following table).
- \Rightarrow Use the **SET** key to confirm the selected unit

	Display	Conversion factor
		1 g =
Gram	g	1.
Ounce	ΟZ	0.035273962
Troy ounce	ozt	0.032150747
Pound	lb	0.0022046226
Tael Hongkong	tlh	0.02671725
Tael Taiwan	tlt	0.0266666
Grain (dependent on model)	gn	15.43235835
Pennyweight	dwt	0.643014931
(dependent on model)		
Momme	(mom)	0.2667
Tola	tol	0.0857333381
Carat (dependent on model)	ct	5
Freely selectable factor *)	FFA	XX.XX

*) Input conversion factor

- As specified above, press repeatedly the MODE key until "FFA" appears.
- ➡ To enter the factor, press the SET key; the enabled digit starts flashing.

Using the **MODE** button, the displayed value is increased by 1, with the **PRINT** button it is reduced by 1. With the **TARE** key selection of the number to the left.

- \Rightarrow Confirm input by pressing the **SET** key.
- ⇒ Press repeatedly the SET key to take over the "Freely selectable factor" as current weighing unit.

Dosing and The Auto-Zero function is used to tare small variations in weight automatically.

In the event that small quantities are removed or added to the material to be weighed, incorrect weighing results can be displayed due to the "stability compensation" in the balance. (Example: Slowly draining fluids from a container on the balance). When apportioning involves small variations of weight, it is advisable to switch off this function.


If **Zero-Tracking** however is switched off, the weighing display becomes more busy.



⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.



- \Rightarrow Press the **MODE** button several times until **"tr"** is displayed.
- - \Rightarrow Acknowledge using **SET** key, the current setting is displayed.

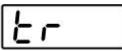

 \Rightarrow Select the desired settings by pressing the **MODE** key.

tr	on	Function activated
tr	off	Function deactivated

- \Rightarrow Use the **SET** key to confirm selection.
- Selection adjustment weight

In the model series KERN FKB_A, the adjustment weight can be selected from three pre-set nominal values (approx.1/3; 2/3; max) (refer also to following table, factory setting with grey background). In order to achieve high-quality weighing results in the sense of the measuring technology, it is recommended to select the nominal value as high as possible. The non delivered adjustment weights can be purchased from KERN as option.



nur



- ⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.
- \Rightarrow Press the **MODE** key several times until **"CAL**" is displayed.
- \Rightarrow Acknowledge using **SET** key, the current setting is displayed.
- \Rightarrow Select the desired settings by pressing the **MODE** key.
 - \Rightarrow Use the **SET** key to confirm selection.

PCB 100-3	CB 100-3 PCB 200-2		PCB 350-3	PCB 1000-1		
20g	50g	50g	100g	200g		
50g	50g 100g		150g	500g		
70g	150g	150g	200g	700g		
100 a	200 a	200g	300g	1000~		
100g	200g	250g	350g	1000g		

PCB 1000-2	B 1000-2 PCB 2000-1		PCB 3500-2	PCB 6000-0	
200g	500g	500g	1000g	1000g	
500g	500g 1000g		1500g	2000g	
700g	1500g	1500g	2000g	5000g	
1000a	2000a	2000g	3000g	6000 a	
1000g	2000g	2500g	3500g	6000g	

PCB 6000-1	PCB 10000-1
1000g	2000 g
2000g	5000g
5000g	7000g
6000g	10000g

Display background illumination



- ⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.
- ⇒ Press the **MODE** button several times until **"bl**" is displayed.
- \Rightarrow Acknowledge using **SET** key, the current setting is displayed.
- \Rightarrow Select the desired settings by pressing the **MODE** key.

bl	on	Background illumination switched on	Contrastful display which can also be red in the darkness.					
bl	off	Background illumination switched off	Battery saving					
bl	Ch	The background illumination will be switched off automatically 10 sec after having reached a stable weighing value.	Battery saving					

 \Rightarrow Use the **SET** key to confirm selection.

Animal weighing function

The animal weighing function can be applied for busy weighing. During a defined period the average value of the weighing results is formed.

The more unquiet the weighed item, the longer the period should be selected.



- ⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.
- ⇒ Press the **MODE** button several times until **"ANL**" is displayed.
- ⇒ Acknowledge using **SET** key, the current setting is displayed.
- ⇒ Select the desired settings by pressing the **MODE** key

ANL	3	Period 3 seconds				
ANL	ANL 5 Period 5 seconds					
ANL	ANL 10 Period 10 seconds					
ANL	ANL 15 Period 15 seconds					
ANL	ANL off Animal weighing not active					

- \Rightarrow Use the SET key to confirm selection.
- Put the weighing good (animal) on the weighing plate and press the SET button. In the display runs a "Countdown". The average value of the weighing results is displayed and remains displayed on the screen.
- ⇒ Use the SET key to change between animal weighing and normal weighing.
- ⇒ Press the SET key to restart the weighing cycle for animal weighing.

Reset to factory setting

This function resets all balance settings to factory setting.



- ⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.
- \Rightarrow Press the **MODE** button several times until "**rSt**" is displayed.
- \Rightarrow Acknowledge using **SET** key, the current setting is displayed.
- ⇒ Select the desired settings by pressing the **MODE** key

rSt	yes	Balance will be reset to factory setting.
rSt	no	The balance keeps its individual setting

⇒ Use the SET key to confirm selection. The balance returns to weighing mode.

9.4 Interface parameters

Data output is carried out via interface RS 232 C.

General

The previous condition for the data transfer between balance and a peripherical device (e.g. printer, PC ...) is that the appliances are set to the same interface parameters (e.g. baud rate, transfer mode ...).

Data transfer mode



- ⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.
- \Rightarrow Press the **MODE** button several times until **"Pr**" is displayed.
- \Rightarrow Acknowledge using **SET** key, the current setting is displayed.
- ⇒ Select the desired settings by pressing the **MODE** key

rE CR	Data output via remote control commands						
Pr PC Data output using the PRINT key							
AU PC	AU PC Continuous data output						
bA Pr	Output on bar code printer						
AU Pr	Autom. data output of stable weighing values						

⇒ Use the SET key to confirm selection. The balance returns to weighing mode.

printout Using this function data are selected which are to be sent via the RS232C (**not** valid for data transfer mode BAPr).



- ⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.
- \Rightarrow Press the **MODE** key several times until **"LAPr**" is displayed.
- \Rightarrow Acknowledge using **SET** key, the current setting is displayed.
- ⇒ Select the desired output parameter by pressing the **MODE** key

Hdr	Edition of the headlines
GrS	Edition of the total weight
Net	Edition of the net weight
tAr	Edition of the tare weight
N7E	Edition of the stored weight
PCS	Edition of quantity
AUJ	Edition of the unit weight
Rqt	Edition of the reference quantity
FFd	Edition of a page feeding at start printer output
FFE	Edition of a page feeding at end printer output

- After actuating the SET button, the current state is displayed (on / off).
- ⇒ Use MODE and PRINT key to change the status "on ≒ off".
- ⇒ Use the SET key to confirm selection. The balance returns to weighing mode.

By that way the user can configurate his own data block, which then is sent to a printer or to a PC.

1

Baud rate

The baud rate defines the transfer speed vie the interface, 1 Baud = 1 Bit/second.



- ⇒ In weighing mode keep the **PRINT** key pressed until **[Unit]** appears.
- ⇒ Press the **MODE** key several times until **"bAUd**" is displayed.
- ⇒ Acknowledge using **SET** key, the current setting is displayed.
- ⇒ Use **MODE** key select the desired settings

9600 ⇒ 4800 ⇒ 2400 ⇒ 1200 ⇒ 19200

⇒ Use the SET key to confirm selection. The balance returns to weighing mode.


10 Data output RS 232 C

10.1 Technical Data

- 8-bit ASCII Code
- 1 start bit, 8 data bits, 1 stop bit, no parity bit
- Baud rate selectable at 1200, 2400, 4800, 9600 and 19200 baud
- Miniature plug-in necessary (9 pole D-Sub)
- For operation with interface faultless operation is only ensured with the correct KERN interface cable (max. 2m)

10.2 Pin allocation of balance output bushing:

Front view:

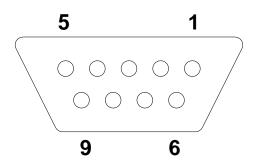

- Pin 2: Transmit data
- Pin 3: Receive data
- Pin 5: Signal ground

10.3 Explanation of the data transfer Pr PC:

Press the **PRINT** key, at stable weight the format is transferred from **LAPR**.

a.	Form	at fo	r stat	ole va	alues	for v	veigh	t/qua	ntity/	/perc	entag	je	

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
М	S	N ₁	N_2	N_3	N_4	N_5	N_6	N_7	N_8	N ₉	N ₁₀	В	U_1	U_2	U_3	CR	LF

b. Format in case of fault

N .			0000	01.10													
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
В	В	В	В	В	В	В	В	В	В	В	E	r	r	0	r	CR	LF

AU Pr:

As soon as the weighing value is stable, the format is automatically transferred from **LAPR**.

c. Format for stable values for weight/quantity/percentage

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Μ	S	N_1	N_2	N ₃	N_4	N_5	N_6	N_7	N_8	N ₉	N ₁₀	В	U_1	U_2	U ₃	CR	LF

d. Format in case of fault

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
В	В	В	В	В	В	В	В	В	В	В	Е	r	r	0	r	CR	LF

AU PC:

The weighing values are sent automatically and continuously, no matter if the value is stable or unstable.

_							<u></u>					-					
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Μ	S	N ₁	N_2	N_3	N_4	N_5	N_6	N ₇	N_8	N ₉	N ₁₀	В	U_1	U_2	U_3	CR	LF

f. Format in case of fault

			0000	01.10													
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
В	В	В	В	В	В	В	В	В	В	В	Е	r	r	0	r	CR	LF

g. Format for unstable values for weight/quantity/percentage

								3									
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Μ	S	N ₁	N_2	N ₃	N_4	N_5	N_6	N ₇	N_8	N ₉	N ₁₀	В	В	В	В	CR	LF

rE Cr:

The remote control commands s/w/t are sent from the remote control unit to the balance as ASCII code. After the balance having received the s/w/t commands, it will send the following data.

Take into account that the following remote control commands must be sent without a subsequent CR LF.

- **s** Function: Stable weighing value for the weight is sent via the RS232 interface
- w Function: Weighing value for the weight (stable or unstable) is sent via the RS232 interface
- t Function: No data are sent, the balance carries out the tare function.

h. Format for stable values for weight/quantity/percentage

1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Μ	S	N_1	N_2	N_3	N_4	N_5	N_6	N_7	N ₈	N ₉	N ₁₀	В	U_1	U_2	U_3	CR	LF

I. FUIMALIII CASE ULIAUIL	i.	Format in case of fault	
---------------------------	----	-------------------------	--

			0000		~~~~												
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
В	В	В	В	В	В	В	В	В	В	В	Е	r	r	0	r	CR	LF

j.	Forn	nat fo	r uns	stable	e valu	ies fo	or wei	ight/c	quant	ity/pe	ercent	tage					
1	2	3	4	5	6	7	8	9	10	11	12	13	14	15	16	17	18
Μ	S	N ₁	N ₂	N ₃	N₄	N ₅	N ₆	N ₇	N ₈	N۹	N ₁₀	В	В	В	В	CR	LF

Symbols

М	Blank or M
S	Blank or minus sign (-)
N ₁ N ₁₀	10 numeric ASCII codes for weight values including decimal places or blanks
$U_1 \dots U_3$	3 ASCII codes for weighing unit pcs. / % / or blank
В	Blank
E, o, r	ASCII code or "E, o, r"
CR	Carriage Return
LF	Line Feed

10.4 Output on bar code printer

The data transfer mode has to be set on **"BA Pr"** (chapter 9.4).

As bar code printer the Zebra printer model LP2824 is provided.

Take into account that the output format of the balance is fixedly defined and cannot be changed.

The printer format is stored in the printer, i.e. in case of a failure the printer cannot be changed with a new one from factory, previously it is necessary that KERN installs the respective software.

The Zebra printer and the balance must be connected to the delivered interface cable when they are switched off.

After switching-on both appliances, and after reaching the status ready-for-operation, a label will be printed out when pressing the **PRINT** button.

11 Service, maintenance, disposal

11.1 CLEANING

Before cleaning, disconnect the appliance from the operating voltage.

Please do not use aggressive cleaning agents (solvents or similar agents), but a cloth dampened with mild soap suds. Take care that the device is not penetrated by fluids and polish it with a dry soft cloth.

Loose residue sample/powder can be removed carefully with a brush or manual vacuum cleaner.

Spilled weighing goods must be removed immediately.

11.2 Service, maintenance

The appliance may only be opened by trained service technicians who are authorized by KERN.

Before opening, disconnect from power supply.

11.3 Disposal

Disposal of packaging and appliance must be carried out by operator according to valid national or regional law of the location where the appliance is used.

12 Instant help

In case of an error in the program process, briefly turn off the balance and disconnect from power supply. The weighing process must then be restarted from the beginning.

Help:

Fault

Possible cause

- The displayed weight does not glow.
- The balance is not switched on.
- The mains supply connection has been • interrupted (mains cable not plugged in/faulty).
- Power supply interrupted.
- Batteries are inserted incorrectly or empty
- No batteries inserted.

The displayed weight is permanently • Draught/air movement changing

- Table/floor vibrations
- The weighing plate is in contact with foreign matter.
- Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

The weighing value is obviously wrong

- The display of the balance is not at zero
- Adjustment is no longer correct.
- Great fluctuations in temperature.
- Electromagnetic fields / static charging (choose different location/switch off interfering device if possible)

Should other error messages occur, switch balance off and then on again. If the error message remains inform manufacturer.