


FR-CTN 60g x 0.001g

Before Weighing

Turn the scale on and leave to warm up for 15 minutes before using, to allow the scale to function correctly and produce accurate results.

Error Messages

EEEE: Overload

: Low battery

Weighing Procedures

1. Press [ON/OFF] to turn on the scale.

When the scale is turned on it does a warm up sequence before showing 0.000g on the display.

2. Selecting a weighing unit

Press [UNIT] to select the unit of measurement you require. You can choose from, g, ct, oz, dwt, ozt, gn, & mg.

3. Start weighing

Without a container

Make sure the scale reads 0 before weighing.

If it does not press the [ZERO] button and wait for the stable indicator to appear on the display (it looks like an eye).

When the stable indicator appears you are ready to begin weighing.

Place the object on the middle of the weighing pan and wait for the stable indicator to appear again before recording the reading.

Weighing with a container

Place your empty container on to the weighing pan and press the [ZERO] button to set the display to 0.

Wait for the stable indicator to appear, and then place your object/s in the container.

Wait for the stable indicator to appear again before recording the reading.

FUNCTION SETTINGS

(To change auto power off and back light settings)

To enter function setting mode, turn the scale off using the [ON/OFF] button.

Press and hold [ZERO] button and keep [ZERO] pressed whilst turning the scale on with the [ON/OFF] button and release the [ZERO] button when either A_ON or A_OFF appears on the display.

Press the [ZERO] button to change A-OFF to A-ON or vice versa.

A-OFF stands for disable auto power off.

A-ON stands for enable auto power off.

Now press [UNIT] to change the display to show L-ON or L-OFF.

Press [ZERO] to change from L-OFF or L-ON or vice versa.

L-OFF stands for disable back light function.

L-ON stands for enable back light function.

Press [ON/OFF] button when you are ready to exit the function settings.

SCALIX
MULTIPURPOSE SCALES

UK Distributor

ATP Instrumentation Ltd

Tournament Way

Ivanhoe Industrial Estate

Ashby De La Zouch

Leicestershire

LE65 2UU

Tel 01530 566800

Fax01530 560373

sales@atp-instruments.co.uk

www.atp-instrumentation.co.uk